



SOP-BCR-4.16

Preparation of PBS

Author: S. Clouthier

Approved: M. Wicha

Rev:
1.0

Issued: 09/24/98
Revised: 7/24/12

1.0 Purpose

The purpose of SOP 4.16 is to prepare sterile PBS.

2.0 Scope

SOP 4.16 is intended to cover all resources, personnel and equipment in the BCR laboratory.

4.0 Materials and Reagents:

No.	Name	Description	Concentration	Storage Location
1.0	PBS	Gibco Phosphate buffered saline	10X	

5.0 Procedure:

- 5.1 Add 2, 500 mL bottles of 10X PBS to carboy.
- 5.2 Then, add 10 L total of distilled water.
- 5.3 Final concentration should be 1X.

6.0 Reference(s)

Not Applicable.

7.0 Procedural Change Description

Revision	Date	Reference	Description of Change
1.0	7/24/12	CL	Updated procedure